



Controller Event Codes and ASCII Commands

<u>Printer Code</u>	<u>Meaning</u>
F	Filling
LL	Low Level Alarm
LH	High Level Alarm
AL	Temp A Low Alarm
AH	Temp A High Alarm
BL	Temp B Low Alarm
BH	Temp B High Alarm
PF	Power Failure Alarm
HG	Hot Gas Bypass Alarm
BV	Low Battery Alarm
US	Usage Alarm
FT	Fill Time Alarm
LO	Lid Open
CA	Temp A Calibration Alarm
CB	Temp B Calibration Alarm
CG	Bypass Temp Calibration Alarm
BY	Bypassing
FD	Filling Disabled
ZO	Level Zeroing
AM	Alarm Muted
BB	Running on battery backup

ASCII Commands

ALMS	INITEE	HITBS?	LOTB?	FILL?	HILM?	RATE?
ALMS?	LNSATP	HITBM	LOTBS?	FILLM	LEVL?	DATE
BAUD	LNSATP?	HITBM?	LOTBM	FILLM?	LFIL	DATE?
BAUD?	BPTMP?	LOGPER	LOTBM?	FILT	LFIL?	TIME
CALTA	HITA	LOGPER?	TEMPA?	FILT?	LOLA	TIME?
CALTB	HITA?	LOTA	TEMPB?	HFIL	LOLA?	UNID
CALVL	HITAS?	LOTA?	TUNI	HFIL?	LOLS?	UNID?
CALVL?	HITAM	LOTAS?	TUNI?	HILA	LOLM	CLEVLG
*IDN?	HITAM?	LOTAM	FILAS?	HILA?	LOLM?	EVNCT?
HISTA	HITB	LOTAM?	FILL 1	HILS?	LUNI	EVENT?
HISTB	HITB?	LOTB	FILL 0	HILM	LUNI?	EVNLOG?

See Chart Connect 3000 Manual for detailed descriptions and usage

Chart Industries, Inc.
 MVE, BioMedical Group
 2200 Industrial Drive, Suite 500
 Ball Ground, GA 30107
 Ph 770-721-7759 • Toll Free 1-800-482-2473
 Fax 770-721-7758



www.chartbiomed.com

Chart BioMedical, Ltd.
 Unit 2, Maxdata Centre
 Downmill Road, Bracknell
 Berks RG12 1QS, United Kingdom
 Ph +44(0) 1344 403100
 Fax +44(0) 1344 429224

Copyright © 2009 Chart Industries