

Having trouble viewing this email? [Click here](#)



## MVE Chart Tech Tips

In This Issue

[Product Information](#)  
[Helpful Hints / FAQs](#)  
[Accessories](#)

PDF Version

To view Tech Tips in PDF,  
click here.

Join Our List


[Join Our Mailing List!](#)

August, 2011

### PRODUCT INFORMATION

#### **MVE Freezer Inventory Systems** Square Racks / Vial Boxes / Blood Bag Frames / Blood Bag Cassettes

Chart MVE freezers are sold without any racks or frames. These components are sold separately. Please contact your distributor for an MVE Freezer rack catalog. The catalog will list part numbers and rack system sizes specifically for Chart freezers. For more information, please go to the [Chart Biomedical Web site](#).

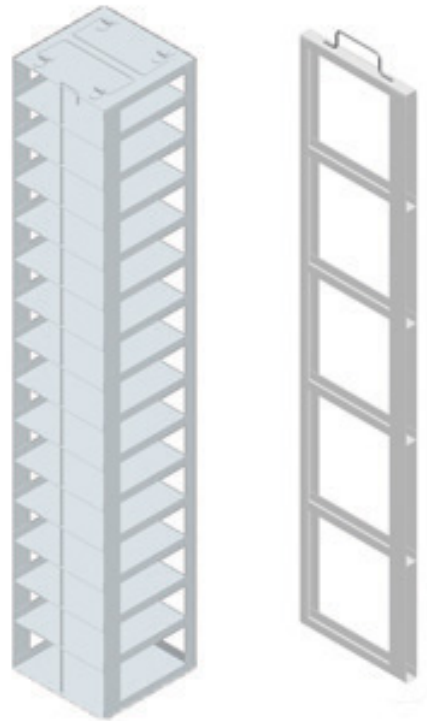

MVE Freezer Racks

#### **MVE Freezer Plumbing Systems**

While performing routine freezer maintenance, be sure to regularly inspect all plumbing and piping connections to ensure no leaks exist,

especially for the copper and brass fittings. Transfer hoses that are frequently removed may potentially leak if not properly secured.



MVE Plumbing Assembly



HE Plumbing Assembly

## MVE Series

The upper chamber temperature of the MVE Series open top freezers is  $-125^{\circ}\text{C}$  ( $-257^{\circ}\text{F}$ ) when used as vapor storage.



MVE Series

## HELPFUL HINTS / FAQs

### Freezer Neck & Lid Assemblies

**Q:** Why are the freezer neck and lid constantly forming ice?

**A:** Extreme humidity levels may cause increased condensation and frost to occur on the freezer neck and lid assemblies. A practical approach and a good prevention method is to wipe down the freezer neck and lid assembly during the maintenance inspection. Inspect the lid gasket for wear and replace as necessary.

**Q:** What is the best method to order a lid gasket?


**A:** Provide the serial and model number to the freezer. At times it may become necessary to provide a photograph and dimensions to ensure accurate replacement of freezer lids.

## ACCESSORIES

### Vial Boxes

Chart sells two 2" and 3" plastic and cardboard boxes for vial storage.

#### Plastic Boxes:

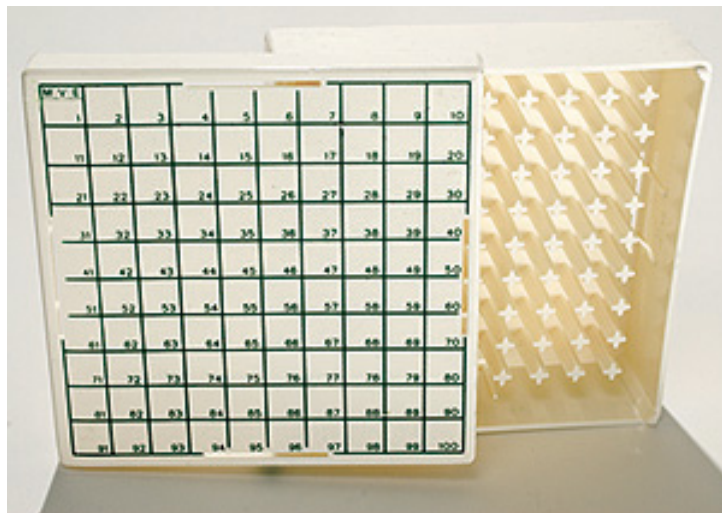

100 Vial Plastic Box

9719649 Holds 100 interior threaded vials 2" high

9719659 Holds 100 interior threaded vials 3" high

#### Cardboard Boxes:





100 Vial Cardboard Box

9714099 Used for 81 external threaded vial-2" high  
9714109 Used for 81 external threaded vial-3" high

## Contact Us

### For ordering information contact Customer Service:

United Kingdom	+44 (0) 1344 403 100
France	+33 (0) 9707 54 646
Germany	+49 (0) 202 739 55429
Italy	+39 049 784 0142
	customerservice.europe@chart-ind.com
Australia/New Zealand	+61 2 9749 4333
Asia/Pac Rim	770.721.7759
The Americas	770.721.7759
US Toll Free	800.482.2473
	customerservice.usa@chart-ind.com

### For technical information contact Technical Service:

United Kingdom	+44 (0) 1344 403 100
France	+33 (0) 5610 82 977
Germany	+49 (0) 202 739 55407
Italy	+39 049 644 498
	techservice.europe@chart-ind.com
Asia/Pac Rim	770.721.7759
	techservice.usa@chart-ind.com
The Americas	770.721.7759
	techservice.usa@chart-ind.com
US Toll Free	800.482.2473
	techservice.usa@chart-ind.com

[Forward email](#)



Try it FREE today.

This email was sent to snystrom@sequal.com by [inga.dolezar@chartindustries.com](mailto:inga.dolezar@chartindustries.com) | [Update Profile/Email Address](#) | Instant removal with [SafeUnsubscribe™](#) | [Privacy Policy](#).

Chart Industries | 2200 Airport Industrial Dr. | Ste. 500 | Ball Ground | GA | 30107