


Fog Clear Lid Switch Harness

The lid switch functionality can be altered for open top freezers. The Fog Clear Lid Switch Harness bypasses the lid switch, allowing the unit to automatically perform a fog clear when the lid is lifted. If a fill cycle is initiated, the freezer will fill while the lid is open. Another nice feature of the fog clear harness is every time the lid is opened, it is documented in the data event log of the controller. Installing the apparatus is straightforward; it simply plugs into the terminating end of the number 3 and 6 pins labeled Lid Switch. The TEC 3000 settings alter the capabilities of the fog clear harness. In all cases, under the add-on menus, the lid switch must be enabled. If Lid Open Fog Clear is disabled, the hot gas bypass time restarts and the fill time continues from the previous count, every time the lid is opened and closed. If Lid Open Fog Clear is enabled on the controller, the hot gas bypass and fill cycle does not stop, regardless if the lid is open or closed. Below, you will find a schematic. If interested, order using part number 13296488 FOG CLEAR LID SW HARNESS.

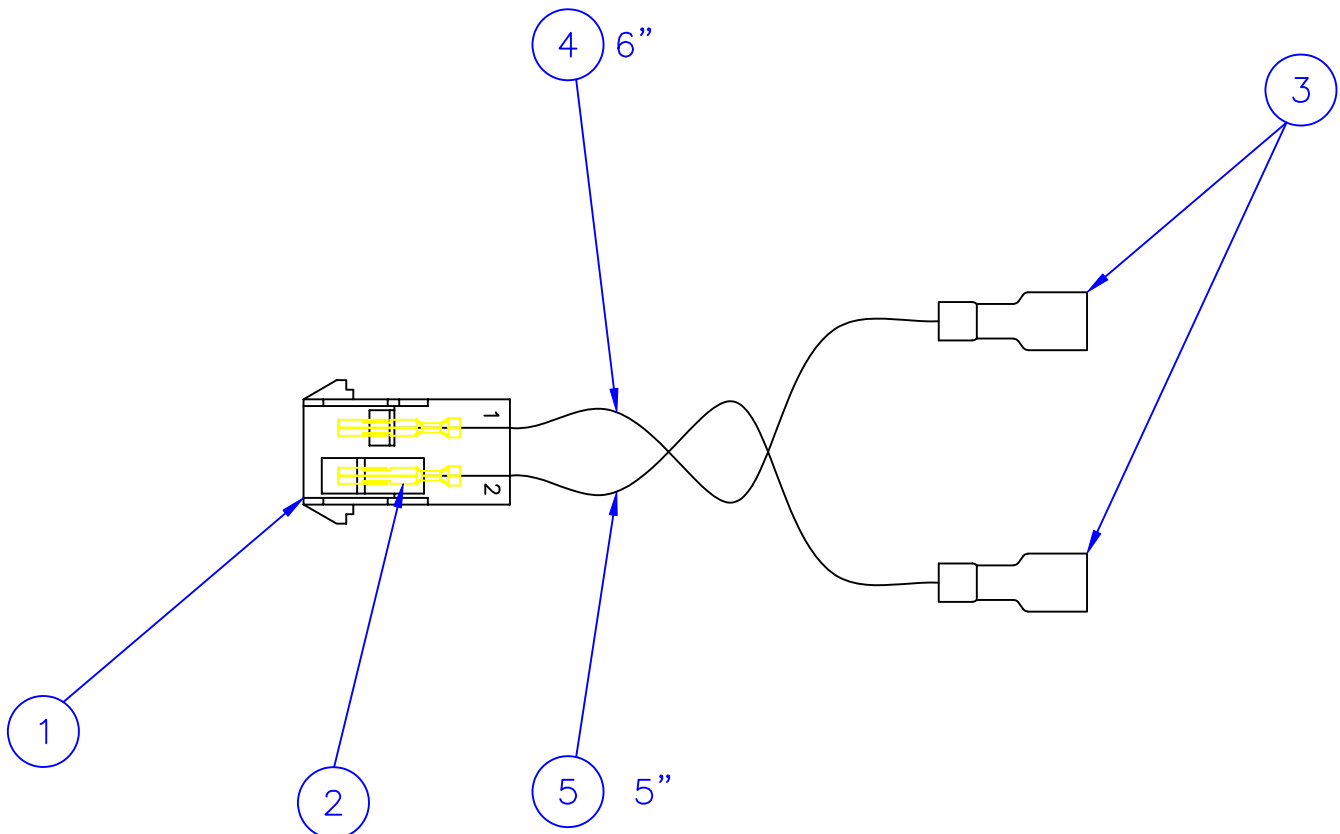

New Face at Chart BioMedical

Please welcome our latest employee to come on board Chart, Holly Hankins. Holly hails from Birmingham, Alabama and enjoys the adventures of being a mom. She will be using her knowledge and experience to serve the BioMedical Division as a Domestic Customer Service Representative. Please make sure to give Holly a warm welcome.



Customer/Technical Info

For ordering information contact

Customer Service:

USA & Canada: 800-482-2473
Europe: +44-1344-403100
Asia/Australia: +61-297-494333

For technical information contact

Technical Service:

USA & Canada: 800-482-2473
LaDahvia Flournoy: ladahvia.flournoy@chart-ind.com
Ramone Gonzalez: ramon.gonzalez@chart-ind.com
Johnny Greene: jonathan.greene@chart-ind.com
UK & Europe: +44-1344-403100

w w w . c h a r t b i o m e d . c o m