



# MVE Tech Tips



November & December, 2008

## Printer Kit Settings

In order to successfully connect a TEC 3000 unit to a printer using the printer kit, settings on both the controller and the printer need to be defined. The controller needs to be in Printer mode and the baud rate set to 9600. The printer should be set on the following

Print Direction .....	Bi-D
Software.....	ESC/P
High speed draft.....	On
I/F mode.....	Serial
Auto I/F wait time .....	10 seconds
Baud rate .....	9600BPS
Parity .....	None
Data Length .....	8bit
Parallel I/F bidirectional mone .....	On
Packet Mode .....	Auto
Character Table .....	PC 437
International character set for Italic table .....	Italic U.S.A.
Auto CR (IBM 2380 Plus).....	Off
IBM character table .....	Table1

The default settings can be changed from the printer's control panel. Make sure paper is loaded into the printer and turn the unit off. Wait 5 seconds, while depressing the Tear Off button, turn on the printer. Pressing the Tear Off button again prints the instructions to change the default settings.

## Cabosil Inserts for a Cryoshipper Mini

---

Below is some information for the cabosil inserts designed to be used in conjunction with the Cryoshipper Mini. The inserts can be mixed and matched to the customers specifications.

Insert Size	Amt of Cabosil (lbs)	Part Number	Hold Time (days)
2"	0.75	11046952	2
3"	1.00	11046944	2.8
4"	1.5	11046936	4

Note: The hold time was determined using 2 inserts

## Technical Tips Distribution List

---

If you know someone who would benefit from receiving the monthly Tech Tips, please send their name, company, and email address to [elizabeth.hofstetter@chart-ind.com](mailto:elizabeth.hofstetter@chart-ind.com).

Please send an email to the same address if you wish to discontinue your monthly subscription to our Technical Tips.

## Customer/Technical Info

---

### For ordering information contact

#### Customer Service:

USA & Canada: 800-482-2473  
Europe: +44-1344-403100  
Asia/Australia: +61-297-494333

### For technical information contact

#### Technical Service:

USA & Canada: 800-482-2473  
LaDahvia Flournoy: [ladahvia.flournoy@chart-ind.com](mailto:ladahvia.flournoy@chart-ind.com)  
Ramone Gonzalez: [ramon.gonzalez@chart-ind.com](mailto:ramon.gonzalez@chart-ind.com)  
Johnny Greene: [jonathan.greene@chart-ind.com](mailto:jonathan.greene@chart-ind.com)  
UK & Europe: +44-1344-403100  
Gary Call: [gary.call@chart-ind.com](mailto:gary.call@chart-ind.com)

**www.chartbiomed.com**